

THE ULTIMATE CHALLENGE:

High-temperature filter for depyrogenation tunnels & ovens.

By Michael Feldtmann



HOW INDUSTRY TRENDS CAN LEAD TO NEW CONCEPTS AND IDEAS FOR HIGH TEMPERATURE FILTRATION.

In the life science industry, hot air up to 350 °C is utilized in static ovens and depyrogenation tunnels to sterilize glass vials, ampoules, and cartridges before they are aseptically filled. The hot air for this process is typically filtered through special high-temperature HEPA filters (HT filters) before it is distributed in a unidirectional airflow through the tunnel.

To avoid product contamination, cleanroom quality air is needed throughout the tunnel. These air filters are therefore of the highest particle collecting efficiency and HT filters of H13 grade are typically used today to filter the air for the hot zone (ISO 5 in the working area acc. to ISO 14644). The choice of the right HT filters is critical to ensure consistent and optimum product safety, as well as maximum production quality and uptime.


The same applies to OEM suppliers of depyrogenation tunnels, who also need effective air filters to guarantee the reliable function and performance of their specialized equipment.

STERILIZATION AND DEPYROGENATION PROCESSES

Manufacturers in the life science industry, such as pharmaceutical companies, use both sterilization and depyrogenation techniques to ensure product quality and safety. Each process has a specific purpose:

- Sterilization renders products free of microorganisms, including bacterial endospores. Sterilization eliminates all forms of life and biological agents. Common methods are the use of high temperatures or intense radiation.
- Depyrogenation removes or inactivates bacterial endotoxin based on a “D-value” – the time required to kill 90% of the exposed microorganisms at a given temperature. The D-value is also used to assess microbial thermal resistance and analyze the thermal death time.

ENDOTOXINS

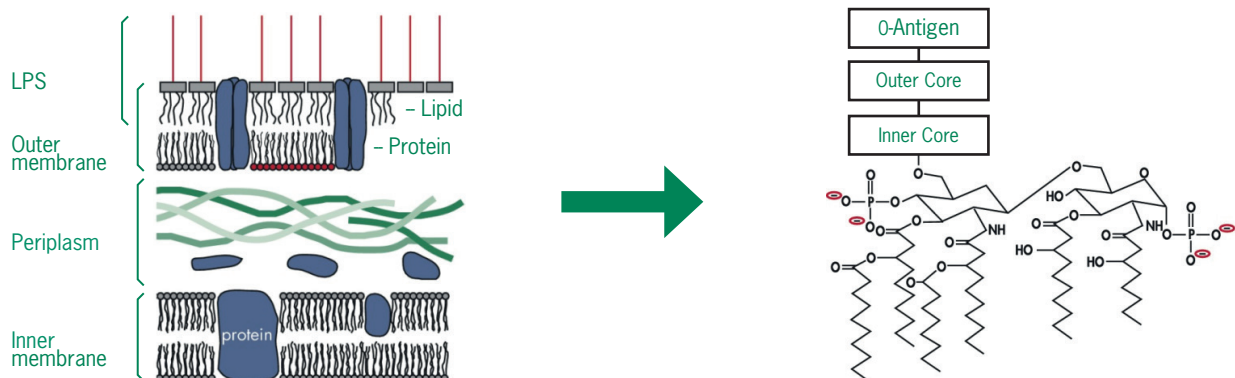

INFLUENCE OF ENDOTOXINS ON BIOLOGICAL APPLICATIONS

Endotoxins strongly influence transfection of DNA into primary cells and sensitive cultured cells, and increased endotoxin levels lead to sharply reduced transfection efficiencies. Furthermore, it is extremely important to use endotoxin-free plasmid DNA for gene therapy applications, since endotoxins cause fever, endotoxic shock syndrome, and activation of the complement cascade in animals and humans. Endotoxins also interfere with in vitro transfection into immune cells such as macrophages and B cells by causing nonspecific activation of immune responses. These responses include the induced synthesis of immune mediators such as IL-1 and prostaglandin. It is important to make sure that plastic ware, media, sera, and plasmid DNA are free of LPS contamination to avoid misinterpretation of experimental results.

WHY HEPA FILTRATION IS NEEDED

In tunnel sterilization systems and depyrogenation tunnels, three important system components have to function optimally to satisfy safety and hygiene standards: the heater elements, the conveyor belt speed and belt recorder, and the HEPA filters. The last are highly important because they ensure cleanroom-quality air and eliminate particles that can contaminate the product and impact production yields.

In the depyrogenation tunnels used in pharmaceutical processing, these high temperature filters are installed directly over the conveyor belt where vials, ampoules or cartridges are transported. The HT filters are specially made to protect these ultraclean processes and have to meet the most stringent requirements and standards. They are designed to maintain integrity and rated performance values at extremely high temperatures.

PROBLEMS ASSOCIATED WITH HT FILTERS

The offering on the market are a wide selection of HT filters for depyrogenation tunnels, but certain filter products do not perform optimally for the application. Some HT filters will release particulate contamination during the high temperature aseptic filling process itself.

This impacts output and quality negatively and results in expensive downtime and filter changes. The particle emission source is often the sealant or other materials used in the filter's construction, which can mechanically interact during periods of rapid temperature change.


INDUSTRY DEMANDS

HT filter customers have expressed four critical needs for cleanroom quality air:

- Ensure the maintenance of ISO Class 5 even when temperature is changing
- Extend HT filter product life time
- Reduce the HT filter installation, preparation, and cleaning time to eliminate the cleaning & preset afterwards and reduce the preparation time for production
- Reduce process downtime.

End users' ultimate goal is to maximize the production yield of processes creating their highest value products. With that in mind the trend is going towards new filter solutions with highly reduced emissions and with no requirement for tempering or cleaning. Future performance requirements for such filter are to meet the highest compliance standards for providing a consistent Grade A manufacturing environment and boosts process yields for tunnels used in the life science industry. The ISO 5 requirement in the operating area is to be the key and further on this the design innovations have to be questioned for filter pack, media and construction. A filter to be installed and integrated into processes up to 350°C without any tempering or cleaning. Reducing start-up and operating costs for the depyrogenation tunnel is key.

In addition, any pharmaceutical-grade filter should have the same efficiency as the other H14-grade filtration in depyrogenation tunnel hot zones, exceeding former HT filter limitations of H13 efficiency.

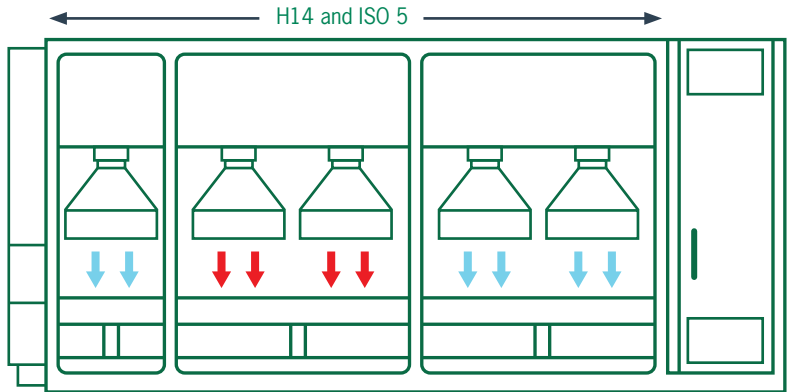

TESTING FOR INTEGRITY AND PERFORMANCE


Due to the crucial function and role of these filters, all filters are factory tested prior to delivery and installation. The measurements conducts leak-tests on every high-temperature filter in a highly controlled clean environment, documenting all test results in accordance with the ISO 9001-2000 quality system.

Enhancing the sampling capability upstream of the filter scanning systems with special dilution equipment is mandatory to measure these high filtration efficiencies. The entire filter face, including the media to frame interface, should be scanned by a specially shaped isokinetic probe traversing in an overlapping pattern and at an appropriate speed to uncover the presence of any leaks. Each Filter should carry an individual scan test report certifying 100% leak free performance. The individual scan test report which is supplied with all Filter are basis for a secured process on customer base.

Di-Ethyl-Hexyl-Sebacate (DEHS), an aerosol liquid used to challenge cleanroom filters during scanning, is released upstream of the filters in a particle size of 0.12 to 0.17 microns to test unit leakage.

The efficiency of HT filters is determined according to EN 1882, the European standard for determining the efficiency of Efficiency Particulate Air (EPA), High Efficiency Particulate Air (HEPA) and Ultra Low Penetration Air (ULPA) filters. This standard determines the ability of the filters to collect and remove the Most Penetrating Particle Size (MPPS).

When comparing the particulate removal capabilities of the H13 and H14 filters the benefits of the higher filter class are obvious. (See Figure X the 10X increase in particle efficiency at MPPS resulting from the substitution of an H14 filter for an H13 unit).

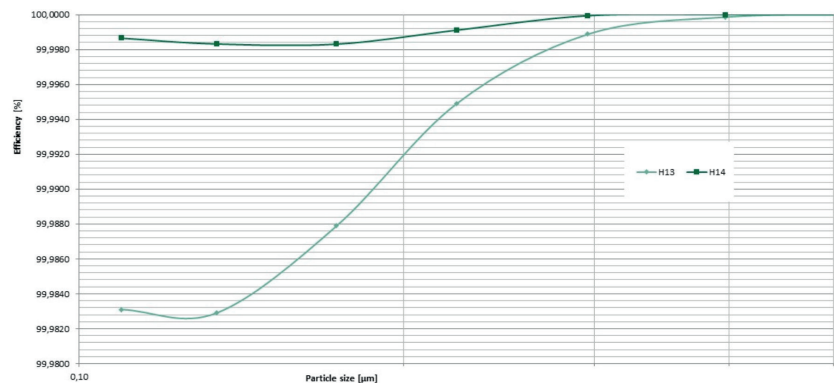

The present depyrogenation tunnel market needs reliable HT filters of H14 efficiency for safety and security of high-temperature production processes. ISO Class 5 cleanroom conditions are required throughout the depyrogenation tunnel. It is therefore extremely important that the HT filters do not emit particles during tempering and temperature changes. They must also remain leak-free during normal operation at steady temperature.

A future design should match benefits including:

- Quality performance throughout all production phases
- To achieve a fast production set up the filter need not to be tempered or prebaked before use, so particulate emissions is reduced.
- To achieve maintenance of < ISO 5 in hot and cool zones of the tunnel, even when temperatures vary, ultra-low emissions should be enabled.
- 350°C working temperature; (Short-term peak temperatures to 400°C.)
- To achieve an extended life time a sealed, airtight filter pack in a reinforced frame could be a solution.
- Filter media free of bisphenol A, phthalates and formaldehyde.
- A reliable H14 filter quality also for the Hot Zone .
- High reliability: leak-free, consistent performance after a large number of cycles (>150-200 cycles).
- With the endurance of temperature fluxes up to +5°C per minute the process requirements of ISO class 5 to be maintained.

An important requirement is No “pre-baking” before production. The filter should be ready to operate immediately after installation and emits no fumes, eliminating the need to maintain an elevated temperature during the suspension of production or when the depyrogenation tunnel is not in operation, saving energy.

New development show excellent results when using for example a resin having a geopolymer type reaction (PSDS) This sealant has been specially formulated to be consistent with the other components of the filter. These polymers have similar properties as inorganic polymers

This geopolymer has properties similar to plastics, but contains no dangerous solvents, does not burn, and does not outgas toxic fumes. Like a refractory, it is resistant to chemical attack and erosion. This type of reaction allows the geopolymer resin to withstand a temperatures up to 600°C

This specially designed, patented geopolymer matrix has a dimensional stability for temperatures exceeding 450°C and will not shrink, or leak. No protection grid is needed.

A new construction should include the newest material developments, specifically stainless steel materials to reduce elongation at higher temperature, ensuring dimensional stability. The stainless-steel separators for the deep pleats also minimize the risk of particle emission during depyrogenation and the sterilization processes.

CERTIFIED PERFORMANCE AND MATERIALS

Such applications require strict compliance with safety, traceability and control. Tested and developed according to precise specifications, filters for all filtration stages should be manufactured with materials that will not contaminate delicate production processes.

Reliability for manufacturers that cannot afford to take risks. The products should be specially designed for life science industries requiring filters to be resistant to decontamination agents, and for food and beverage industries needing components certified for food contact. The materials in the filters also have to be inert against microbial growth, but at the same time, they must be free of any harmful chemical components.

All filters should be compliant with FDA, REACH, EC1935:2004, VDI6022 and ISO 846 raw-material specifications. Letters of conformance should accompany every filter to meet product compliance requirements.



ABOUT THE AUTHOR

Michael Feldtmann Wirtschaftsing. BA

E-mail: michael.feldtmann@camfil.com

Web: www.camfil.com

Michael Feldtmann is the Product Manager for Camfil, the world leader in air filtration solutions and filter production. His responsibility covers the Product Management of the Clean Process in Europe, which includes all airfilter products with very high Efficiency and the belonging housing systems. He is working in the air filtration business for over 30 years, responsible for different product portfolios. His profession is Industrial Engineer, Dual Study in cooperation with the Drägerwerk AG in Lübeck. As the german expert in the ISO TC 142 WG 4 (HEPA and ULPA Filter) he is belonging to the german and the international standardisation process for High Airfiltration purposes. As part of his job he has worked over several years with all types of High Temperature Hepa filter and the needs from the industry.

Camfil – a global leader in air filters and clean air solutions

For more than half a century, Camfil has been helping people breathe cleaner air. As a leading manufacturer of premium clean air solutions, we provide commercial and industrial systems for air filtration and air pollution control that improve worker and equipment productivity, minimize energy use, and benefit human health and the environment.

We firmly believe that the best solutions for our customers are the best solutions for our planet, too. That's why every step of the way – from design to delivery and across the product life cycle – we consider the impact of what we do on people and on the world around us. Through a fresh approach to problem-solving, innovative design, precise process control and a strong customer focus we aim to conserve more, use less and find better ways – so we can all breathe easier.

The Camfil Group is headquartered in Stockholm, Sweden, and has 30 manufacturing sites, six R&D centres, local sales offices in 30 countries, and 4,500 employees and growing. We proudly serve and support customers in a wide variety of industries and in communities across the world. To discover how Camfil can help you to protect people, processes and the environment.

www.camfil.com

